



385PR

**G-Biosciences ♦ 1-800-628-7730 ♦ 1-314-991-6034 ♦ [technical@GBiosciences.com](mailto:technical@GBiosciences.com)**

A Geno Technology, Inc. (USA) brand name

# Prediluted Gamma-Globulin Standard

(Cat. # 786-114G)



think proteins! think G-Biosciences [www.GBiosciences.com](http://www.GBiosciences.com)

## INTRODUCTION

Prediluted Gamma-Globulin protein standards are in solution made in saline buffer containing 0.05% sodium azide. They are provided in six different concentrations ranging from 0.1mg/ml to 1.0mg/ml.

## ITEM(S) SUPPLIED (Cat. # 786-114G)

Description	Size
Prediluted Gamma Globulin 0.1mg/ml	5ml
Prediluted Gamma Globulin 0.2mg/ml	5ml
Prediluted Gamma Globulin 0.3mg/ml	5ml
Prediluted Gamma Globulin 0.5mg/ml	5ml
Prediluted Gamma Globulin 0.8mg/ml	5ml
Prediluted Gamma Globulin 1.0mg/ml	5ml

## STORAGE CONDITION


The kit is shipped at ambient temperature. Store refrigerated at 4°C, upon arrival. For long term storage keep at -20°C and is stable for 1 year. Avoid frequent freeze-thaw cycle.

## PROTOCOL

For use with protein estimation kits, use 50µl from each vial for 1.5ml assay format and 5µl per well for 96-well plate assay. For specific types of protein assays, use an appropriate volume of the standard as needed.

## RELATED PRODUCTS

Download our Protein Assays Handbook.

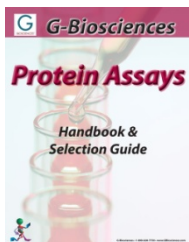

<http://info.gbiosciences.com/complete-protein-assay-guide>

For other related products, visit our website at [www.GBiosciences.com](http://www.GBiosciences.com) or contact us.

*This page is intentionally left blank*



**[www.GBiosciences.com](http://www.GBiosciences.com)**