

VWR® PCR PLATE SPINNER

Unique

Compact

Economical




VWR® PCR PLATE SPINNER



- Space saving design
- Easy plate loading system
- Quiet, simple operation
- Cold-room compatible

The VWR PCR Plate Spinner offers a personal sized centrifuge designed to spin down PCR plates. The unique centrifuge is equipped with a dual position vertical rotor that accepts a wide variety of PCR plates; 96-well, 384-well, skirted and non-skirted plates are all compatible with the Plate Spinner. The rotor features integral plate guides, ensuring that the plates stay vertical and do not tip inside the rotor. With the use of the available adapter plate (89184-612) you can spin down individual PCR tubes, 8 or 12 place PCR strips, and up to 4 x 48 well half plates.

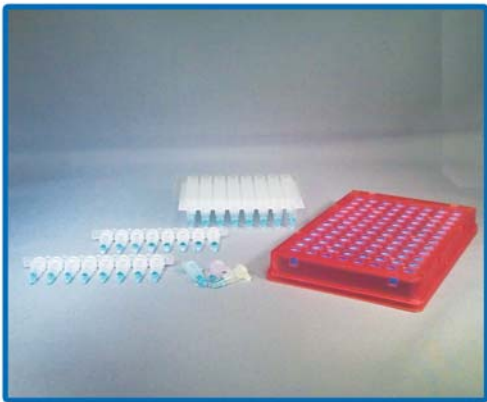

Sealed microplates are loaded vertically through the loading slots. The rotor can be turned using the rotor knob to insert the second plate. Upon closing the lid of the unit, the motor quickly accelerates to 2,500 RPM and will maintain speed until the OPEN button is pressed. After pressing the button, the rotor brake engages and brings the rotor to a quick gentle stop.



Specifications

Maximum Speed	2,500 RPM
Maximum RCF	500 x g
Capacity	2 PCR Plates
Dimensions (W x D x H)	19 x 21 x 18.3 cm 7.5 x 8.2 x 7.2 in
Weight	3 kg / 6 lbs
Electrical	120V or 230V, 50/60 Hz

Description	Cat No.
VWR PCR Plate Spinner, 120V/60Hz, with 2 Plate Rotor	89184-608
VWR PCR Plate Spinner, 230V/50z, with 2 Plate Rotor	89184-610
Adapter for 8 x 0.2 mL PCR Strips, 12 x 0.2 mL PCR Strips, 0.2 mL PCR Tubes, or 4 x 6 Half Plates	89184-612



1.800.932.5000 | vwr.com

Prices and product details are current when published; subject to change without notice. | Certain products may be limited by federal, state, provincial, or local regulations. | VWR makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC. All prices are in US dollars unless otherwise noted. Offers valid in US and Canada, void where prohibited by law or company policy, while supplies last. | VWR, the VWR logo and variations on the foregoing are registered (®) or unregistered trademarks and service marks, of VWR International, LLC and its related companies. All other marks referenced are registered by their respective owner(s). | Visit vwr.com to view our privacy policy, trademark owners and additional disclaimers. ©2016 VWR International, LLC. All rights reserved.