

TMB Microwell Peroxidase Substrate System



| <u>Catalog No.</u> | <u>Size</u> |
|--------------------|-------------|
| 50-76-00 | 600 mL |
| 50-76-03 | 2700 mL |
| 50-76-11 | 200 mL |

DESCRIPTION

TMB Microwell Peroxidase Substrate System is a 2 component system which develops a deep blue color when reacted with horseradish peroxidase conjugates in ELISA. It is not recommended for blotting or immunohistochemical assays. See KPL's catalog for appropriate substrates.

FORM

Catalog No. 50-76-00 consists of the following:
3 x 100 mL TMB Peroxidase Substrate (Catalog No. 50-76-01)
3 x 100 mL Peroxidase Substrate Solution B (Catalog No. 50-65-00)
Catalog No. 50-76-03 consists of the following:
3 x 450 mL TMB Peroxidase Substrate (Catalog No. 50-76-02)
3 x 450 mL Peroxidase Substrate Solution B (Catalog No. 50-65-02)
Catalog No. 50-76-11 consists of the following:
1 x 100 mL TMB Peroxidase Substrate (Catalog No. 50-76-01)
1 x 100 mL Peroxidase Substrate Solution B (Catalog No. 50-65-00)

STORAGE/STABILITY

Store both components at 2 - 8°C. Stable for a minimum of 1 year from date of receipt when stored at 2 - 8°C. The TMB Peroxidase Substrate solution may develop a yellow tinge over time which does not affect product performance.

CONTENT

TMB Peroxidase Substrate contains 3,3',5,5'- tetramethylbenzidine at a concentration of 0.4 g/L in an organic base. Peroxidase Substrate Solution B contains H₂O₂ at a concentration of 0.02% in a Citric Acid buffer.

USE

Preparation:

Mix equal volumes of TMB Peroxidase Substrate and Peroxidase Substrate Solution B in a clean, preferably HDPE, polypropylene or glass container immediately prior to use. Substrate solution should remain clear. Warm to room temperature before use.

Substrate Development:

Following incubation with peroxidase labeled conjugate, wash plate thoroughly. Add 100 µL prepared substrate solution to each well. As the color develops, tap gently to mix. Incubation times will vary depending on your assay.

To Stop Reaction:

Stop reaction by adding an equal volume of stop solution to the microwell plate. This will halt color development and will turn the TMB substrate yellow. (See Recommended Stop Solution).

To Read Reaction:

After stopping, read at a wavelength of 450 nm. Stopped reaction should be read within 30 minutes.

When to Stop Substrate Reaction:

Upon addition of stop solution, absorbance values increase 2 - 3 fold. The point at which the substrate reaction is stopped is often determined by the ELISA reader. The O.D. values of the plate should be monitored and the reaction stopped before positive wells are no longer recordable.

To Reduce Substrate Intensity:

High background and/or precipitate in the wells are a sign of overreaction with TMB. To reduce the intensity of the substrate reaction, dilution of the primary antibody and/or conjugate is recommended. Dilution of the substrate is not recommended.

ABSORBANCE MEASUREMENTS

Kinetic Assays: The TMB substrate produces a blue color upon reaction with peroxidase. Read at a wavelength of 620 - 650 nm.

Endpoint Assays: The addition of 100 µL (or an equal volume) of stop solution to the microwell plate will halt color development and will turn the TMB substrate yellow. Read at a wavelength of 450 nm. Stopped reactions should be read within 30 minutes.

RECOMMENDED STOP SOLUTION

The addition of an equal volume (100 µL) of 1M Phosphoric Acid (H₃PO₄) provides the most stable endpoint color. Other commonly used acids (HCl, H₂SO₄) are less efficient in stopping color development.

NOTE: The addition of TMB Membrane Enhancer, Catalog No. 50-77-01, allows the TMB Microwell Substrate to be used on membranes.

PRODUCT SAFETY AND HANDLING

See MSDS (Material Safety Data Sheet) for this product.

TMB Microwell Peroxidase Substrate System

| <u>Catalog No.</u> | <u>Size</u> |
|--------------------|-------------|
| 50-76-00 | 600 mL |
| 50-76-03 | 2700 mL |
| 50-76-11 | 200 mL |

RELATED PRODUCTS

| | |
|--|----------------------|
| Milk Diluent/Blocking Solution Concentrate | Catalog No. 50-82-01 |
| 10% BSA Diluent/Blocking Solution | Catalog No. 50-61-00 |
| Detector [™] Block | Catalog No. 71-83-00 |

The product listed herein is for research use only and is not intended for use in human or clinical diagnosis. Nothing disclosed herein is to be construed as a recommendation to use this product in violation of any patents. The information presented above is believed to be accurate. However, said information and product are offered without warranty or guarantee since the ultimate conditions of use and the variability of the materials treated are beyond our control. We cannot be responsible for patent infringements or other violations that may occur with the use of this product. No claims beyond replacement of unacceptable material or refund of purchase price shall be allowed. All claims regarding product performance must be made within 30 days following date of delivery.

Limited Use License

The purchase of this product conveys to the buyer the non-transferable right to use the product in research conducted by the buyer. The buyer cannot sell or otherwise transfer this product or materials made by use of this product to a third party or otherwise use this product or materials made with this product for Commercial Purposes without written approval of KPL, Inc. For additional information, please visit www.kpl.com or the KPL catalog. To obtain a license or approval to use this product for purposes other than those permitted above, contact Director of Sales at (301) 948-7755.